



## Implementation of HTRF<sup>®</sup> Assay Technology on Infinite<sup>®</sup> F200 PRO

Reader Control Kit, cAMP HiRange Kit, TNF-alpha Kit

### Abstract

The present application note summarizes the results of experiments with the HTRF (Homogeneous Time-Resolved Fluorescence, Cisbio Bioassays) technology on Infinite F200 PRO.

### Introduction

HTRF combines time-resolved fluorescence (TRF) and fluorescence resonance energy transfer (FRET). It is based on the energy transfer between two fluorophores, a long-lived europium or terbium cryptate donor and either a chemically modified allophycocyanine (XL665) or a d2 acceptor. HTRF-based applications have become increasingly popular for the analysis of various molecular interactions and binding studies because of their homogeneity, robustness, sensitivity, and potential for miniaturization.

Tecan's Infinite F200 PRO is equipped with a sophisticated and ingenious mirror system containing a dichroic mirror with a break point at 510nm, which permits optimal excitation in the UV range and thus enables HTRF measurements in the Infinite F200 PRO.

#### HTRF Reader Control Kit

The HTRF Reader Control Kit (RCK) is intended for the calibration of HTRF compatible instruments and for validation of their ability to perform HTRF measurements. The RCK contains the complete analytical chemistry that is necessary to assess the HTRF performance of the reader to be tested and can be used to compare different instruments.

#### cAMP HiRange Kit

Cyclic AMP (cyclic adenosine 3', 5'-monophosphate) is one of the most important intracellular messengers and is involved in numerous biological processes such as GPCRs, protein kinase activation or ion channel regulation. The cAMP HiRange kit is designed for the direct quantification of cAMP in suspension or adherent cells by means of a competitive immunoassay on the basis of HTRF technology. It permits the measurement of agonist and antagonist effects on GPCRs in different cell lines. The measured TR-FRET signal is inversely proportional to the cAMP concentration in the sample.

### TNF-alpha Kit

Tumor necrosis factor (TNF)-alpha is a cytokine involved in systemic inflammation and is a member of a group of cytokines that stimulate the acute phase immune reaction. Under inflammatory conditions TNF-alpha is secreted by activated macrophages and monocytes. The Human TNF-alpha kit allows for the quantitation of TNF-alpha in cells or supernatants by antibody-mediated recognition of two distinct epitopes that are tagged with HTRF donor and acceptor fluorophores. The HTRF ratio increases linearly with the TNF-alpha concentration.

## Material and methods

### Instrument

- Infinite F200 PRO filter-based detection system, equipped with a 510 dichroic mirror

### Microplates

- 96-well half area microplates, white (Greiner Bio-One, Germany)
- 384-well small volume microplates, white (Greiner Bio-One, Germany)

### Reagents

- HTRF Reader Control Kit (62RCLPEA, Cisbio)
- cAMP HiRange kit (62AM6PEB, Cisbio)
- Human TNF-alpha kit (62TNFPEB, Cisbio)

### Assay protocol

#### HTRF Reader Control Kit

The HTRF RCK was used according to the manufacturer's instructions. The measurement was performed in a white 96-well half area plate using 100 µl filling volume per well. The plate was measured in triplicates.

	1	2	3	4	5	6	7	8	9	10	11	12	
A	620 nm Control	Standard 0	Low Control	High Control	Buffer Blank	XL665 Blank	Cryp Blank						
B													
C													
D													
E													
F													
G													
H													

Figure 1: Plate layout for the HTRF Reader Control Kit

- (A1-D1): 50 µl 620 nm control, 50 µl reconstitution buffer
- (A2-H3): 50 µl diluent, 25 µl cryp, 25 µl XL665
- (A4-H4): 50 µl low cal., 25 µl cryp, 25 µl XL665
- (A5-H5): 50 µl high cal., 25 µl cryp, 25 µl XL665
- (A6-C6): 50 µl diluent, 50 µl reconst. buffer
- (A7-C7): 50 µl diluent, 25 µl XL665, 25 µl reconst. buffer
- (A8-C8): 50 µl diluent, 25 µl cryp, 25 µl reconst. buffer

RCK - Passed/failed criteria

Signal-to-Blank ratio	≥ 40 = passed
CV Standard 0	≤ 10 % = passed
CV high Standard	≤ 10 % = passed
CV low Standard	≤ 10 % = passed
Delta F low	≥ 15 = passed
Delta F high	≥ 600 = passed

Table 1: Performance criteria for HTRF certification

### cAMP HiRange Kit

The cAMP HiRange kit was used according to the manufacturer's instructions. The measurement was performed in a white 384-well small volume plate using 20 µl filling volume per well. The plate was measured in triplicates.

	1	2	3	4	...	...	...	...	...	...	...	24
A	Standard 8	Cryp blank										
B	Standard 7											
C	Standard 6											
D	Standard 5											
E	Standard 4											
F	Standard 3											
G	Standard 2											
H	Standard 1											
...												
P												

Figure 2: Plate layout for the cAMP HiRange Kit

- (A1-A3): Standard 8 = 2800 nM cAMP
- (B1-B3): Standard 7 = 700 nM cAMP
- (C1-C3): Standard 6 = 174 nM cAMP
- (D1-D3): Standard 5 = 43.75 nM cAMP
- (E1-E3): Standard 4 = 10.94 nM cAMP
- (F1-F3): Standard 3 = 2.73 nM cAMP
- (G1-G3): Standard 2 = 0.86 nM cAMP
- (H1-H3): Standard 1 = 0.17 nM cAMP
- (A4-C4): Cryptate Bank = cryptate conjugate buffer

### Human TNF-alpha Kit

The Human TNF-alpha kit was used according to the manufacturer's instructions. The measurement was performed in a white 384-well small volume plate using 20 µl filling volume per well. The plate was measured in triplicates.

	1	2	3	4	...	...	...	...	...	...	...	24
A	Calibrator 7											
B	Calibrator 6											
C	Calibrator 5											
D	Calibrator 4											
E	Calibrator 3											
F	Calibrator 2											
G	Calibrator 1											
H	Calibrator 0											
...												
P												

Figure 3: Plate layout for the Human TNF-alpha Kit

**(A1-A24):** Calibrator 7 = 2000 pg/ml TNF-alpha

**(B1-B24):** Calibrator 6 = 1000 pg/ml TNF-alpha

**(C1-C24):** Calibrator 5 = 500 pg/ml TNF-alpha

**(D1-D24):** Calibrator 4 = 200 pg/ml TNF-alpha

**(E1-E24):** Calibrator 3 = 100 pg/ml TNF-alpha

**(F1-F24):** Calibrator 2 = 50 pg/ml TNF-alpha

**(G1-G24):** Calibrator 1 = 20 pg/ml TNF-alpha

**(H1-H24):** Calibrator 0 = diluent

### Measurement settings

Measurement 1 (Donor)	
EX wavelength	320 (25) nm
EM wavelength	620 (10) nm
Mirror	dichroic 510
Lag time	150 µs
Integration time	500 µs
Number of flashes	50
Gain	optimal
Measurement 2 (Acceptor)	
EX wavelength	320 (25) nm
EM wavelength	665 (8) nm
Mirror	dichroic 510
Lag time	150 µs
Integration time	500 µs
Number of flashes	50
Gain	optimal

Table 2: Optimized instruments settings for HTRF measurements on Infinite F200 PRO

### Calculations

#### Detection limit (TNF-alpha):

Detection limit = dose of mean zero + 2 SD

#### Delta R:

Delta R = Average Standard X – Average Std 0

#### Delta F:

$$\text{Delta F} = \frac{\text{Delta R} * 100}{\text{Average Std 0}}$$

#### Assay window (cAMP):

$$\text{Assay window} = \frac{\text{Delta F}_{\text{Std 0}}}{\text{Delta F}_{\text{Std 8}}}$$

## Results

### HTRF Reader Control Kit

The measurement results obtained with the Reader Control Kit demonstrate the applicability of the Infinite F200 PRO for HTRF analyses. The Infinite F200 PRO meets all performance criteria specified by Cisbio (tables 1 and 3), resulting in an exceptionally good measurement range (Delta F low = 32, Delta F high = 1189, Z' value = 0.95) and minimal measurement variations (CVs of all calibrators and standards ≤ 2%).

### Reader Control Kit – Measurement results

Measurement time	2 min 59 s
Gain 620 / 665	162 / 182
Signal-to-Blank ratio	268
CV Standard 0	1.24 %
CV Low Calibrator	1.26 %
CV High Calibrator	1.35 %
Delta R low	278
Delta R high	10454
Delta F low	32
Delta F high	1189
Z' value	0.95

Table 3: RCK measurement results

### cAMP HiRange Kit

As shown in figure 4, the Delta F values obtained with the cAMP dilution series are inversely proportional to the cAMP concentration, resulting in the sigmoidal shape of the curve that is typical for competitive assays.

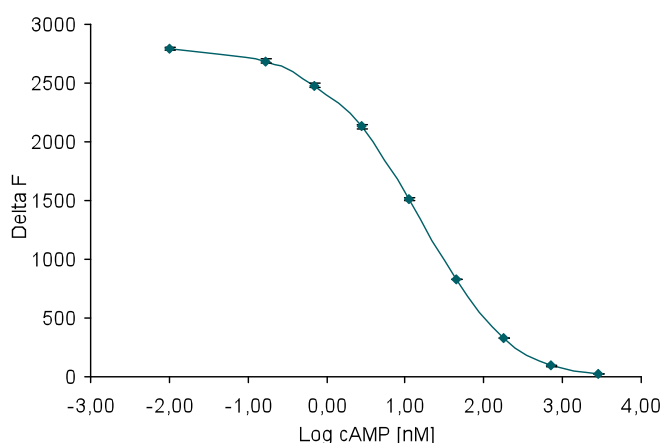


Figure 4: Competitive cAMP curve

### cAMP HiRange Kit – Measurement results

Measurement time	1 min 58 s
Gain 620 / 665	144 / 152
CV Standard o	3.44 %
Delta F Standard 8	20
Delta F Standard 7	94
Delta F Standard 6	330
Delta F Standard 5	827
Delta F Standard 4	1514
Delta F Standard 3	2130
Delta F Standard 2	2480
Delta F Standard 1	2689
Delta F Standard o	2794
Assay window	143

Table 4: cAMP measurement results

### Human TNF-alpha Kit

The measurement of a dilution series of the TNF-alpha standard shows a nearly linear course ( $R^2 = 0.9998$ ) from 2000 to 20 pg/ml TNF-alpha (figure 5).

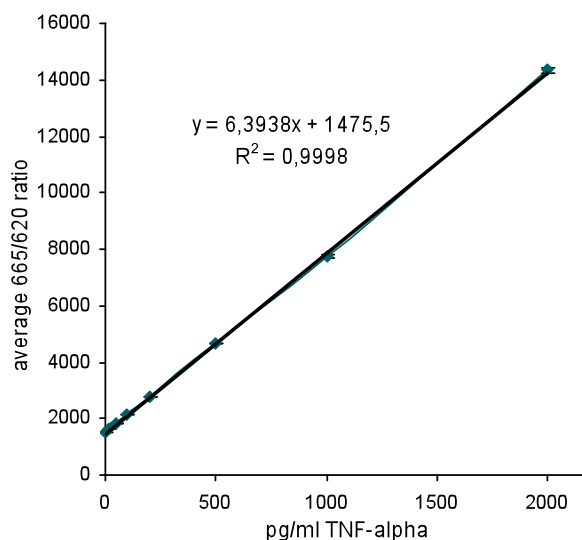


Figure 5: Linearity of TNF-alpha dilution series

According to the evaluation protocol provided by Cisbio, only calibrators 1 to 3 were taken as a basis for the calculation of the TNF-alpha detection limit. The resulting detection limit for TNF-alpha was calculated to be 5,56 pg/ml. The detailed results (average of three independent measurements) are summarized in table 5.

### Human TNF-alpha – Measurement results

Measurement time	9 min 45 s
Gain 620 / 665	162 / 194
CV Standard o	1.26%
Delta F Calibrator 3	41
Delta F Calibrator 2	21
Delta F Calibrator 1	8
Linearity (R2)	0.9998
Detection limit	5,56 pg/ml

Table 5: TNF-alpha measurement results

## Conclusion

In accordance with the criteria for the HTRF Reader Control Kit, the Infinite F200 PRO in combination with the 510 dichroic mirror meets all performance parameters that are relevant for HTRF-compatible readers and is thus regarded as HTRF-certified. In addition, both cAMP and TNF-alpha can be reliably and sensitively quantified in the Infinite F200 PRO.

In summary, the results of the Reader Control Kit, the cAMP HiRange Kit and the Human TNF-alpha Kit show that the Infinite F200 PRO with an integrated 510 dichroic mirror is a perfectly suited detection instrument for the measurement of various HTRF assays formats.

## List of abbreviations

cAMP	cyclic adenosine monophosphate
CV	coefficient of variation
FRET	fluorescence resonance energy transfer
TNF	tumor necrosis factor
TRF	time-resolved fluorescence
TR-FRET	time-resolved fluorescence resonance energy transfer
HTRF	homogenous time-resolved fluorescence
RCK	Reader Control Kit
S:B ratio	signal-to-blank ratio
SD	standard deviation

## Acknowledgements

Tecan would like to express thanks to Marie-Laure Lebreton and Marielle Mazille from Cisbio Bioassays for their valuable comments and support during instrument validation.

.....  
**Austria** +43 62 46 89 33 **Belgium** +32 15 42 13 19 **China** +86 10 5869 5936 **Denmark** +45 70 23 44 50 **France** +33 4 72 76 04 80 **Germany** +49 79 51 94 170  
**Italy** +39 02 92 44 790 **Japan** +81 44 556 73 11 **Netherlands** +31 18 34 48 174 **Portugal** +351 21 000 82 16 **Singapore** +65 644 41 886 **Spain** +34 93 490 01 74  
**Sweden** +46 31 75 44 000 **Switzerland** +41 44 922 89 22 **UK** +44 118 9300 300 **USA** +1 919 361 5200 **Other countries** +41 44 922 8125  
 .....

Tecan and Infinite are registered trademarks of Tecan Group Ltd. Männedorf, Switzerland.

HTRF® is a registered trademark of Cisbio Bioassays. The fluorescence ratio associated with the HTRF® readout is a correction method developed by Cisbio Bioassays that is covered by the US patent 5,527,684 and its foreign equivalents, for which Cisbio Bioassays has granted a license to Tecan Group Ltd. Its application is strictly limited to the use of HTRF reagents and technology, explicitly excluding any other TR-FRET technologies.

© 2010, Tecan Trading AG, Switzerland, all rights reserved.

[www.tecan.com](http://www.tecan.com)